



First-aid plan

Eye injuries

SUBSTANCE	EYEWASH	CONTINUOUS RINSING	RINSING TIME
Foreign bodies Dust, dirt, metal and wood chips etc.	Plum Eye Wash	Plum Eye Wash	Rinse until the foreign body has been rinsed out.
Acids	pH Neutral	Plum Eye Wash	Rinse for approximately two minutes with pH Neutral and continue rinsing with Plum Eye wash until the inconvenience disappears or until you reach a doctor.
Alkali	pH Neutral	Plum Eye Wash	Rinse for approximately two minutes with pH Neutral and continue rinsing with Plum Eye Wash until you reach a doctor.
Other chemicals	Plum Eye Wash or pH Neutral	Plum Eye Wash	Rinse until the inconvenience disappears or until you reach a doctor.
Chemicals on your skin	pH Neutral	Plum Eye Wash	Continue rinsing until the inconvenience disappears.



pH Neutral



Plum Eye Wash



1. Remove the bottle from the station



2. Twist the lid until the sealing breaks.



3. Lean your head backwards and rinse.

Important!

The eye wash fluid must always come in contact with the eye in a gentle, even flow. Press the bottle gently without squeezing it.