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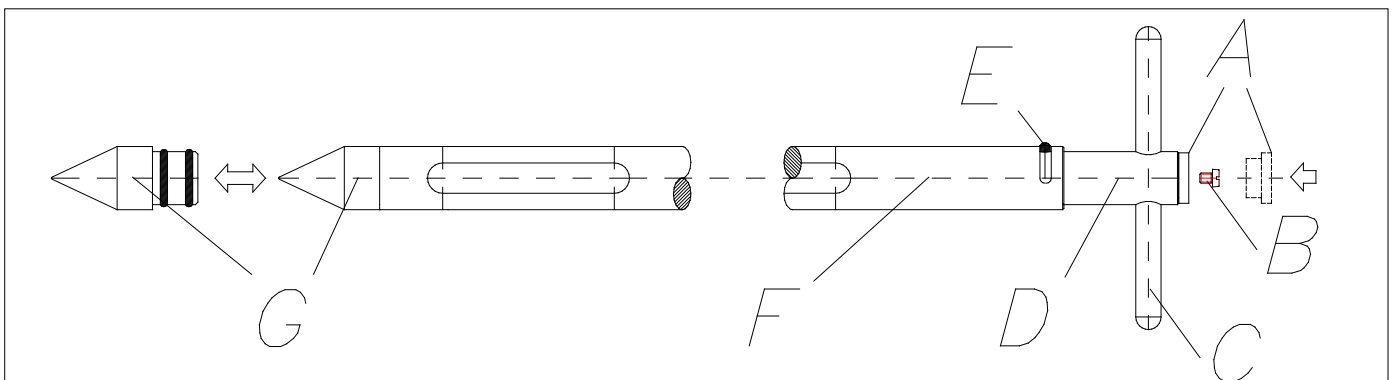
Read carefully before use and keep for later reference

Safety Regulations

Every user must know these instructions and have them at hand at all times. The sampler may only be used for easy-flowing granulates. Do not use the sampler for explosive powders/granulates. This device can only be used within the limits set by the resistance of the materials it is made out of. Sampler parts coming into contact with the sample material are Sampler No. 5317- anodised aluminium/stainless steel Sampler No. 5328- and 5329- stainless steel 1.4305

Follow the reagents suppliers' safety recommendations. Always use the sampler in such a way as not to endanger yourself or others. When using the sampler, never use force.

Packaging/Contents

Please check the contents to see if they are complete or if they have been damaged during transport. Damage to the packaging which can be seen on the outside should be immediately shown to the deliverer/forwarding agent who in turn must confirm it.

Function, Precautionary Measures, Use Restrictions

The device has been designed to take samples of easy-flowing powders or granulates from sacks, big-bags, canisters, barrels. In the case of aggressive substances, protective clothing must be worn, particularly eye and hand protection. The sampler cannot be used for chemicals that attack 1.4305 stainless steel or the anodised aluminum face.

Assembly

Before bringing the sampler into service, the handle should be assembled. For this put the handle "C" through sampling part "D" and fasten the locking screw "B" with a screwdriver. Afterwards introduce the color coding (see chapter "Color Coding").

Color Coding

The coding buttons "A" (red, blue, green, yellow, black) which are supplied together with the sampler, can be used to equip the sampler with a color code, thus preventing that a sampler is accidentally used to sample another substance. The buttons can be easily inserted in the gaps of both handles' ends of the sampling part (see diagram "A"), and are then knocked into their final position using a firm surface.

Sampler Instructions

Set the zone sampler in the "closed" position by rotating the handle clockwise to the right. Insert the zone-sampler to the desired depth in the powder/granulate. Rotate the handle anti-clockwise to the left (now the powder/granulate will flow into the zone-sampler). Now turn the handle clockwise to the right, thus sealing the sample chambers. Carefully remove the zone-sampler and empty the sample onto an appropriately prepared surface.

Warning! Please note that when turning the handle, no parts of the body (finger, palm of the hand) are in the opening. The opening has sharp corners that can cause injury.

Maintenance and Care

The samplers need no maintenance, but particularly after sampling abrasive materials, they should be cleaned as follows:

1. Push in pin "E", grasp handle "C".
2. Pull out "D" completely. To open the outer tube "F", unscrew the tip "G".
3. With the cleaning brush (accessory), the sampling tube can be cleaned thoroughly.
4. Wipe both tubes – inside and out – before re-assembling. The space between inner and outer tube is where jamming can occur.

The samplers should not be lubricated, because this would contaminate the samples. All types of samplers (No. 5317- aluminum/stainless steel or Art.-No. 5328- and 5329- stainless steel models), can be sterilised with steam in the autoclave at temperatures of up to 121 °C.

Accessories/Spare parts

Cleaning brush, ideal zum Reinigen des Sammelrohres 5304-0100

Warning! Avoid injury through chemicals! Wear plastic apron, safety goggles and protective gloves!

Protective clothing

Safety goggles 2502-1001
 Protective gloves 2507-0001
 Plastic apron 2507-1001

Self-adhesive seals for closing your paper sacks

close-it control seal, printed, black 5303-0011
 close-it control seal, printed, red 5303-0012
 close-it control seal, printed, yellow 5303-0014
 close-it control seal, printed, green 5303-0016
 close-it control seal, printed, white 5303-0002

Sample container, sterilisable, made of PP

Screw-cap tubs 100 ml 7103-1000
 Screw-cap tubs 60 ml 6203-0060
 Screw-cap tubs 180 ml 6203-0180

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