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User Manual

Mode d'emploi

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PS 611

PS 620

PS1211

PS1220

P04

P22

P44

DENIOS AG

Dehmer Str. 58-66
32549 Bad Oeynhausen
Tel.: +49 5731 753-0
Fax.: +49 5731 753-199
E-Mail: info@denios.de

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8 General instructions and safety instructions

The instructions in this user manual must be carefully followed and adhered to. All those involved in the installation, operation, maintenance and repair of the product must have read and understood this user manual.

We reserve the right to make technical changes to the illustrations and details in the user manual in order to improve the storage system.

	NOTE
	<ul style="list-style-type: none"> ▪ The user is responsible for installation and commissioning of the acid and alkalis cabinet. ▪ The national guidelines and safety regulations with respect to hazardous materials, safety regulations, occupational safety and operator duties must be observed.

8.1 Intended use of the equipment

- The acid and alkali cabinet is suitable for passive storage indoors.
- Cabinets may only be installed on flat, firm surfaces.

	NOTE
	<ul style="list-style-type: none"> ▪ Do not locate in Ex-zones. ▪ Temperature range -20°C to +40°C. ▪ Do not locate cabinets in direct sunlight.

8.2 Safety instructions

	WARNING
	<ul style="list-style-type: none"> ▪ No fire, naked flames or smoking. ▪ Do not exceed the load capacity of the shelves. ▪ Store materials so that all containers and the spill pallet are visible.

	CAUTION
	<ul style="list-style-type: none"> ▪ There is a danger of crushing when operating the door.

8.2.1 Intended use of the equipment

Acid and alkali cabinets are suitable for the safe, legally compliant storage of water-polluting substances with a flashpoint $>100^{\circ}$ in containers. The volume of an individual container must not exceed the permitted containment volume of the acid and alkali cabinet. The total contents of the stored containers must not exceed ten times the permitted containment volume of the cabinet.

	The storage of flammable liquids with a flashpoint $<100^{\circ}\text{C}$ is not permitted.
--	---

Liquids in DIBt media list 40-1.1 and liquids, mentioned in the following groups of substances, do not require any special proof of resistance:

- Aqueous solutions of organic acids up to 10%
- Mineral acids up to 20 %
- Acidic hydrolysing salts in aqueous solution ($\text{pH}<6$), apart from hydrofluoric acid
- Oxidising acids and salts
- Inorganic alkalis as well as alkali hydrolysing salts in aqueous solution ($\text{pH} > 8$), except for ammonia solutions and oxidising solutions of salts (eg hypochlorite),
- Solutions of non-oxidising salts with a pH value between 6 and 8.

	Containers with water-polluting liquids of various compositions and properties may only be stored together, if it can be proven that these substances will not cause a dangerous reaction.
	When storing media which fall under the hazardous substances regulations, TRGS 510 must be observed.

	NOTE
	<ul style="list-style-type: none"> ▪ The safety instructions in the safety data sheet for the stored substance must be observed.

9 Safety instructions

	WARNING
	<ul style="list-style-type: none"> ▪ No fire, naked flames or smoking. ▪ Do not exceed the load capacity of the shelves. ▪ Store materials so that all containers and the spill pallet are visible.

	CAUTION
	<ul style="list-style-type: none"> ▪ There is a danger of crushing when operating the door.

10 Technical data

10.1 Dimensions and data

Model	W [mm]	D [mm]	H [mm]	Max. load [kg]	Shelves	Spill pallet capacity (l)
PS 611	600	600	1130	50 kg / per shelf	2	35
PS 620	600	600	1990	50 kg / per shelf	4	35
PS 1211	1200	600	1130	50 kg / per shelf	2	75
PS 1220	1200	600	1990	50 kg / per shelf	4	75
P04	460	460	560	20 kg / per shelf	2	5
P22	900	560	870	20 kg / per shelf	2	2x5
P44	900	560	1650	20 kg / per shelf	4	4x5

11 Transport, location and assembly

11.1 Transport

- Only transport the cabinets when empty.

11.2 Set-up conditions

- The acid and alkali cabinets must only be set up on flat, firm surfaces.
- The cabinets must not be set up in a vehicle.
- Ensure the cabinets are protected from damage by vehicles using suitable means (eg impact protection).

11.3 Installation and placement

- The acid and alkali cabinets must be set up level at the final place of use.
- When correctly installed, the wing doors should open and close easily.
- Fit the shelves supplied (depending on design either stainless steel, galvanised steel or plastic) at the required height on the supports in the cabinet.

	CAUTION
	<ul style="list-style-type: none">▪ The PolyStore PS acid and alkali cabinet must be secured to a suitable load bearing wall with the supplied mounting kit. The mounting kit must be adapted as necessary to the actual conditions on site.

11.4 Technical ventilation (accessories)

- Fit the fans in accordance with the instructions supplied separately.

12 Operation

- Unlock the door lock using the key supplied and unbolt the doors by turning the handle to the right through 90°
- Open the cabinet doors
- Check containers are stable after setting them down.
- Close the cabinet by turning the handle back to the horizontal position.
- If necessary, secure the cabinet against unauthorised entry by turning the cylinder lock.

13 Maintenance and Care

- Keep the spill pallet dry and free from contamination.
- Check the spill pallet regularly for leaked fluids.
- Safely dispose of leaked fluids immediately.
- If the spill pallet is damaged, the hazardous materials cabinet must be emptied immediately and taken out of operation.
- Lubricate the locks and hinges with commercially available grease.
- Regular maintenance and care will extend the life of the hazardous materials cabinet.

14 Disposal

	NOTE
	Clean the components thoroughly before disposal to remove any hazardous material residue.

The environmental cabinets and the acid and alkali cabinets are made from various components. After decommissioning, follow local legislation when disposing of any waste.

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